

Post Stand Instruction Manual



Please read the Instruction Manual carefully before installation and keep it for future use.

I. Confirm your purchase order

76MM Post Stand	PZ Post Stand	39MM Post Stand	83MM Post Stand	85MM Post Stand
<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>

II. Open the package and confirm the parts






1	Base
1	Post
1	Post Stand Extension Bar
1	Post Stand Cap
1	Guard Ring
1	Through Hole Focus Rack Check your ordered Through Hole Focus Rack <input type="checkbox"/> 76mm Through Hole Focus Rack <input type="checkbox"/> 83mm Through Hole Focus Rack <input type="checkbox"/> 85mm Through Hole Focus Rack <input type="checkbox"/> 39mm Through Hole Focus Rack <input type="checkbox"/> PZ Microscope Stand
1	Wrench Key
1	95 mm black and white plastic or glass frosted plate (on the Base)
1	Instruction Manual




Open the package and take out the parts carefully. Contact the local dealer for any missing part.

Please check the received parts based on your purchase order.

III. Installation

				
1. Place the Base on the stable table, and thread the post into the base.	2. Lock the side screws with the Wrench Key	3. Thread the Post Extension Bar in the post.	4. Put the Guard Ring on 1/3 of Post Stand, and tighten the handle.	5. Put the Focus Holder on the Post Stand till the Guard Ring and lock the screw knob.

	<p>6. Thread the cap on the post top.</p>		<p>7. Place the black and white plastic or glass frosted plate in the whole by pressing the spring on side</p>
---	---	---	--

Make sure whether the Post Stand you choose includes illumination Base or not. If it excludes illumination Base, please follow the above instruction to install. If it includes illumination Base, please refer to the Instruction Manual and Notes of Illumination Base for using the stand with Illumination Base.

IV. Specifications and Accessories

Post Stand Diameter	1.26" (32mm)
Height of Main Post Stand	12.01" (305mm)
Height of the Post Stand Extension Bar	5.9" (150mm)
Max. Height of Post Stand	17.91" (455mm)
Focusing Working Distance	50mm
Max. Distance from the Post Stand to the front far side of the Base	10.75" (273mm)
Base Dimension	12.6x12.01x0.63" (320x305x16mm)
Can be used with the illumination Base (The Illumination Base is optional.)	

Types of Stands	Types of Microscopes to be matched with the different stands
76MM Post Stand	Can be applied to LEICA, OLYMPUS, NIKON Stereo Microscopes, and etc.
PZ Post Stand	PZ Microscopes
39MM Post Stand	Can be applied to MZ and MZ-II series Microscopes
83MM Post Stand	Can be applied to 83mm Microscopes
85MM Post Stand	Can be applied to MEIJI Microscopes and 85mm Microscopes.

Notes:

1. Save the Wrench Key for your future use. The measurement units for Wrench Key and screws are metric.
2. Please refer to the Instruction Manual and Notes of Illumination Base for using the stand with Illumination Base.
3. Use soft cleaner to clean the metal part of the post, and keep it clean.
4. The working temperature for operation and storage of track stand is from 15°C-45°C.
5. Contact the local dealer for any missing part and problem.