

Track Stand Instruction Manual



Please read the Instruction Manual carefully before installation and keep it for future use.







I. Confirm your purchase order

| | | | | | | |
|--------------------------|--------------------------|--------------------------|--------------------------|--------------------------|--------------------------|--------------------------|
| 76MM Track Stand | PZ Track Stand | 39MM Track Stand | 83MM Track Stand | 85MM Track Stand | B&L Track Stand | N Type Track Stand |
| <input type="checkbox"/> | <input type="checkbox"/> | <input type="checkbox"/> | <input type="checkbox"/> | <input type="checkbox"/> | <input type="checkbox"/> | <input type="checkbox"/> |

Notes: Accessories of the Track Stand with fine-focus are the same.

II. Open the package and confirm the parts

Open the package and take out the parts carefully. Contact the local dealer for any missing part. Please check the received parts based on your purchase order:

| Quantity | Name of Parts | |
|----------|--|--|
| 1 | Base | |
| 4 | Locking screws of the track | |
| 1 | Holder Check your ordered holder |  <input type="checkbox"/> 76mm Holder <input type="checkbox"/> 83mm Holder <input type="checkbox"/> 85mm Holder |
| |  <input type="checkbox"/> 39mm Holder |  <input type="checkbox"/> PZ Microscope Holder |
| |  <input type="checkbox"/> B&L Holder |  <input type="checkbox"/> N Adapter |
| 1 | Track Cap  | |
| 3 | Locking screws of the holder | |
| 1 | Track | |
| 1 | Wrench Key | |
| 1 | 95 mm black and white plastic or glass frosted plate (on the Base) | |
| 1 | Instruction Manual | |

III. Installation

Make sure whether the Track Stand you choose includes illumination Base or not. If it excludes illumination Base, please follow the following instruction to install. If it includes illumination Base, please refer to the Instruction Manual and Notes of Illumination Base for using the stand with Illumination Base.



1. Take out the four screws on the track.



2. Place the Base on an even table by the side, use the Wrench Key to fasten the four screws on the bottom of the track.



3. Take out the three screws on the focus part.



4. Install the microscope Support Holder on the focus setting, and tighten down the three screws.



5. Place the black and white plastic plate in the whole by pressing the spring on side



6. Cover top with cap.

Lay the microscope Stand on an even table after installation.
Put the microscope into the holder, and fasten the screws.

IV. Notes

Notes:

1. Save the Wrench Key for your future use. The measurement units for Wrench Key and screws are metric.
2. Please refer to the Instruction Manual and Notes of Illumination Base for using the stand with Illumination Base.
3. Use soft cleaner to clean the metal part of the post, and keep it clean.
4. The working temperature for operation and storage of track stand is from 15°C-45°C.
5. Contact the local dealer for any missing part and problem.